

# Vernier Bluetooth Radio (BLUE-USB) for use with the Vernier Wireless Dynamics Sensor System

Applies to Ezurio BRBLU03-010A0-02 only



This Bluetooth radio is for use with the Vernier Wireless Dynamics Sensor System (WDSS). It allows communication between the WDSS and a computer running *Logger Pro 3*. No additional interface is required to use the WDSS, although *Logger Pro* will collect data from other devices simultaneously with WDSS collection. Follow instructions below to connect the radio to your computer's USB port. For information on using the WDSS in *Logger Pro*, see the booklet accompanying the WDSS.

No separate installation CD is required.

## System requirements

Windows: Windows XP SP2 or newer<sup>1</sup>, Windows Vista 32 or 64, *Logger Pro 3.7* or newer.  
Macintosh: OS X 10.4, 10.5 or 10.6, *Logger Pro 3.6* or newer.

## Add this Bluetooth radio to your Windows XP or Vista computer

*Logger Pro* requires the Microsoft Bluetooth driver software. The drivers are normally part of Windows XP SP2 and later, but may not be present on your computer if it has not been updated. As a result you may need to connect to the Internet while connecting the radio. You may install *Logger Pro 3.7* or newer either before or after these steps.

1. Connect the PC to the Internet.
2. Connect the Bluetooth radio to an available USB port.
  - a. Windows may display a series of bubbles detailing the installation process. In this case the installation will be complete when the bubbles close.
  - b. Instead of installation bubbles the Windows New Hardware wizard may run. If you are prompted to search the Internet for drivers, accept this search. Allow the Wizard to complete the installation.

## Add this Bluetooth radio to your Macintosh computer

You may install *Logger Pro 3.6* or newer either before or after connecting the radio.

1. Connect the Bluetooth radio to an available USB port.

**See the WDSS manual or visit [www.vernier.com/bthelp](http://www.vernier.com/bthelp) for further information.**

### Vernier Software & Technology

13979 S.W. Millikan Way • Beaverton, OR 97005-2886  
Toll Free (888) 837-6437 • (503) 277-2299 • FAX (503) 277-2440  
[info@vernier.com](mailto:info@vernier.com) • [www.vernier.com](http://www.vernier.com)

Revised 8/2/11



<sup>1</sup> Earlier versions of Windows XP can be updated to SP2 by running Windows Update.